## What's New in iVision-Mac 4.5.0

iVision 4.5 is the first Intel-native version of iVision. It includes the following changes and features:

# **Faster Processing on Intel**

Here are the times for a simple image processing script that reads 20 indexed files, changes them, and writes them back to a different folder:

	G5 (2.3GHz)	MacPro (3GHz)
iVision 4.0	19 seconds	21
iVision 4.5	18	10

Note that camera frame rates are not noticeably improved. The faster processing speeds may speed up scripts that do other processing during the capture loop, though.

# **PCIe support on Intel**

Cameras which use PCI-Express (PCIe) boards are now supported on Intel-based machines.

# Scripts and non-device preferences are not convertible.

Sorry about that.

# File/Image names: 255 Unicode characters

Unicode allows the use of "international" (non-Roman alphabet) characters (e.g. Japanese).

There are some items where Unicode is still not fully supported including:

Row and Column Labels

Drawing Text

To fully support these a new file format will have to be developed, because the older IPLab format does not store these items in Unicode.

# New dialog window name field & resizable dialogs.

The window name field in dialogs is now a single edit field with a menu to the right of it. The menu includes "relative" window names that let you refer to windows based on their position, for use in scripts. The relative names are drawn in italics. To put in a specific name (e.g. "DAPI Image"), pick from the menu, or just type into the field and any italic name will go away.

	Histogram		
Image: Front V	Vindow		
Minimum Bin:	• Use Data Minimum	On:	Front Window Second Window Back Window
Maximum Bin:	• Use Data Maximum	• ROI	DAPI Image FITC Image

Since window names can potentially be very long now, dialogs that contain window name fields can be resized horizontally to see the full name.

	Histog	ram
Image: Project	ABC12345 Study 19254 07-	21-2009 Infusion 4 FITC @ 20x
Minimum Bin:	• Use Data Minimum	On: C Entire POL
Maximum Bin:	• Use Data Maximum	
Number of Bins	(Max - Min) + 1	
		Cancel OK

#### More commands allow you to specify the image to use.

All commands that reference image windows now include a window name field, rather than only working with the front image. This should make scripting easier, since you don't have to worry as much about which image is at the front.

## Open As replaced by enhanced Open command.

The "Open As" command let you tell iVision what format a file was (e.g. TIFF or PICT), so if the file had the wrong extension or file type associated with it you could still open the file. Since it was not used much and seemed to confuse people more than it helped them, the "Open As" has been removed. It's functionality has been included in the "Open" command.

To keep "Open" simple to use, the default file format is set to "Automatic". This will figure out the format of the file for you and open the image if it is in any format that iVision knows.

The "Automatic" format should be used in almost all cases.

The best reason to use a specific format is when using a script to open a large number of images that are all the same format but do not have an extension or other file type (e.g. from an old OS 9 CD). In that case telling iVision the format will speed up the file reading.



## **Open Multiple Files at once, interactively.**

The Open command allows you to select more than one file at a time interactively. This is not scriptable.



## File Sets are just Indexed and Folder (no more List).

#### Segmentation sets gone.

You can just place the segmentation command into the script without creating a segmentation set. This required that the number of colors saved for the Color Cube feature be limited to a maximum of 20.

#### Pseudocolor dialog expanded:

The "Predefined" menu is reduced to the "complex" color ramps, like Rainbow, Classify, and System, as well as the commonly used "Monochrome" and "Reverse Mono".

There are also two new options to create simple color ramps:

Pseudocolor	Pseudocolor
Image: Front Window	Image: Front Window
<ul> <li>Wavelength:</li></ul>	<ul> <li>Wavelength:</li></ul>

The menu to the right of the ramp choices lists "single color" color definitions (e.g. Red, Green, Blue, Cyan, etc.). You can add to this list with the new "Define Colors" command.

## **Edit CLUT replaced with Define Colors:**

The Define Colors command lets you specify a single color and give it a name:

Color:	New	My Orange *Blue	ŕ
As Darkfield Ramp: Apply	Revert Save Save As	*Coomassie Blue *Cyan *Dapi-Bfpschst-Amca *Ethidium Bromide	
As Brightfield Ramp: Apply	Delete	*Fitc-Gfp-Cy2-DiO *Green *Magenta *Red	- and a second
Frame: 🚹 0 📕		*Tritc-Cy3-Dil *TxRed-Pl	

The dialog show what a darkfield and brightfield ramp made from this color will look like. If there is an image open when this dialog is open, you can apply either of the ramps to the image to see how they look. You can also step through the frames of the image to preview what the change will look like.

The colors defined in this command are also available in FCV when assigning colors to images in the Color Join command.

# <u>Changes to</u> Variables

For the most part variables still work the same as before. You still use a number with the "Var" button in a dialog to specify a variable.

But the Variables window has changed from an image window to a palette:

$\Theta \cap \Theta$	Variables
Remove	
Name	Value
v0	6.00
v1	5.00
v3	0.00
v4	42.00
v16	20.00
v32	40.00
v33	1.00
v34	0.00
v200	0.00
v201	0.00
v202	4032.00
v203	2688.00
v204	4032.00
v205	2688.00
v206	4032.00
v207	2688.00
v232	New Window Name
v249	0.20
v255	0.20
v300	6.00
•	

In this window the name of the variable is preceded by a lower case "v".

The Variable window only shows those variables that have been used at some point (and have not been removed). You can change the value of a variable in this window by clicking on the Value column and typing a new value. In addition, variables can contain text as well as numbers.

#### Updated Set Variables dialog.

The Set Variables dialog includes several ways to place text into variables:

- A new "Text" option to set any arbitrary text string.
- A "Window Name" option to place the name of an existing window into a variable.
- The "Acquire Info" menu now includes non-numeric fields such as Filter Names and dates, which will be placed into the variable as a text string.



# Updated Enter Variables dialog.

The Enter Variables dialog has been updated to add some formatting features, and to allow text to be entered into variables:

00	Enter Variables				
Title: Ente	er Values				
Description: Plea	se enter the requested values:				
Usage:	Prompt:		Variable to use:		Display Constant:
Text 🛟	Name of Experiment ?	A V	0		0
Line	]	Å V	1		0
Title 🛟	Camera Settings	A V	2		0
Number 🛟	Exposure?	A Y	10	V	0.200
Number 🛟	Binning?	A Y	11	☑	2
Filter Name 🛟	CRI Filter	];	5		0
✓ Unused		A Y	6		0
Number Text		A Y	7		0
Checkbox		Å	8		0
Title		A V	9		0
			Cano	el	ОК

When the script is run, the result looks like this:

Please enter t	Enter Values he requested values	:
Name of Ex	periment ?	
Contr	ol	
	Camera Settin	gs
Exposure?		0
Binning?		2
CRI Filter:	Red	\$
Script: Untitled :	1	
	C	Cancel OK

#### **Using Variables in Text Fields**

To use any variable (numeric or text) in a text field, you use a special formatting rule that lets you insert the value of the variable into the text.

*Place the name of the variable inside braces:* {*v123*}.

You can optionally use one of two formatting options for numbers: zero fill, or decimal places.

For zero fill you can specify a total length of 2 to 6 digits.

For decimal places you can specify 0 through 6 digits after the decimal place.

You specify the option by following the variable name with a comma and then a 'z' or a 'd', followed by the number.

## **Example:**

if v42 = 12.54700000

[v42}	=> 12	(truncate)
[v42,z2}	=> 12	(truncate)
[v42,z3}	=> 012	(truncate)
[v42,z4}	=> 0012	(truncate)
[v42,z5}	=> 00012	(truncate)
[v42,z6}	=> 000012	(truncate)
[v42,d0}	=> 13	(round)
[v42,d1}	=> 12.5	(round)
[v42,d2}	=> 12.55	(round)
[v42,d3}	=> 12.547	(round)
[v42,d4}	=> 12.5470	(round)
[v42,d5}	=> 12.54700	(round)
[v42,d6}	=> 12.547000	(round)

You can put more than one variable substitution in a text field. For example, if v10 = 4, and v12 = 32.4112:

"Experiment {v10,z3} temperature: {v12,d1} F" would produce: "Experiment 004 temperature: 32.4 F".

Variable substitution is available in:

Window Names	
Draw Text:	text
Enter Variables:	title, description, prompt(s) [for Enter Var and Title uses]
Script Query:	title, message, button titles
Script Alert:	title, message

# Scripts:

The Script window has been updated. It is resizable horizontally.

Record OFF	Co	ntinue Run Q	8
Command		Comment -	
<b>Dispose All Windows</b>			
Show/Hide Variables	-		
Use File Set		Open, var33	
Use File Set		Save, var34, ask	
Delay		start timer	
Set Variable	-	v16 = 0	
Repeat			
Open	-		
Exit If			
Change Color Type	-	up to C48	
Split Color Channels	-	get R, G, B front to back	
Median Filter	-	3x3 on Red	
Change Window	-	Red to back	
Autosegment	-		
Binarize	-		
Change Window	-	Green to back	
Pixel Arithmetic	-		
Change Window	-	Blue to back	
Dispose Window	-	original	
Merge Color Channel	s 🗖		
Rename Window	-	index with v16	
Set Variable	-	v16++	
Save As		to File Set "Results"	
Dispose All Windows			
End Repeat			
Show/Hide Variables	-		
Delay		elapsed seconds into v86	
END			

The END symbol now acts like a label,: you can jump to END without adding a label just above it.

The Comment column can be toggled to show the parameters in the command:



At the moment this is more confusing than helpful because the description shows every possible field even if they aren't in use, and in a hard to read format. This will be improved in the future.

The search field at the top of the window will highlight any commands that contain the given text:



In this example the File Set named "Iris Frames" is used by two commands "Use File Set" and "Open". The command name, the user comment, and the parameters are all searched for the given text. The search field remembers the last 10 searches:



The script commands palette includes two new commands: Enable and Disable.



These commands allow you to disable commands in a script, which causes them to be skipped over, as if they weren't in the script at all. Disabled commands are displayed in red:



## Scripts remember different size/placement than interactive dialogs.

If you are scripting a command where you want the dialog to be opened while the script is running (with the "+"), the size and position of the dialog used when the script is run will be saved independently from the size and position used when not running a script. So you can reliably place the dialog at a specific place on screen.

New File Extensions for iVision files:				
iVision Script		.ivsc		
iVision Preference		.ivpr		
iVision Extension	.ivex			
iVision Device Module	.ivdm			

## **About Box:**

The About Box contains a popup menu which shows the name and version of all Extensions in use. This does not include device modules, which are listed in the Device Setup dialog.

At the bottom of the popup menu is "**Reveal in Finder**", which will open a Finder window for the folder which contains the currently running iVision. This helps verify which copy of iVision is running (if there is more than one on the machine),



The Device Setup dialog has an **Open Folder** button to show the current Device Modules folder.

0	0	0	Device Setup	
	5	Cameras	<u></u>	Setup
	<b>⊠</b>	Hamamatsu DCAM	v 4.5.0 d64	Connect
	R	Roper PVCam	v 4.5.0 d64	Connect
	R	QImaging QCam	v 4.5.0 d64	
	KQ	Scion	v 4.5.0 d64	
	~	Virtual Camera	v 4.5.0 d64 💙	Connect All
	15	Stages		
	$\Delta$	ASI Stages	v 4.5.0 d64	
	Δ	Leica DM6000 Stages	v 4.5.0 d64	
	$\Delta$	Leica Two Stages	v 4.5.0 d64	
	Â	Ludl Stages	v 4.5.0 d64	
	Â	Nikon E1000 Stage	v 4.5.0 d64	
	$\Delta$	Nikon Remote Focus	v 4.5.0 d64	
	$\Delta$	Nikon TE2000 Stage	v 4.5.0 d64	
	$\Delta$	Olympus BX Stage	v 4.5.0 d64	
	$\triangle$	Olympus IX Stage	v 4.5.0 d64 👻	
	Deta	ils:		Tools:
	DB	Ready to Connect.		Ones Falder
	Fai	iled to initialize the Hama	Open Folder	
				Serial Status
	ls t	the camera connected an	v 4.5.0 d64	
				Done
		· · · · · · · · · · · · · · · · · · ·		